

INSTRUCTION LEAFLET**MONOCLONAL Ab POLYPESTI****1ml vial (→ 1:50)**

Mixing of non-conjugated anti-pestivirus monoclonal antibodies for viral revelation tests (BVD-MD, BD and HCV strains).

COMPONENTS

Polyvalent mixing of purified ascitis fluids obtained from mice inoculated with different hybridoma directed against the following pestivirus (glyco-) proteins :
structural envelope gp48, 53, 62, non structural p80-125.

INDICATION OF USE

Immunological revelation of pestiviruses, mainly in tests using the Indirect Immunofluorescence technique (IIF) (i.e : viral isolation).

PROCEDURE

Note : This reagent must be aliquoted (without dilution) and kept deep-frozen in order to ensure an optimal preservation.

Before each use, allow one aliquote to thaw and dilute 1:50 in PBS.

Then use, according to the viral revelation technique processed.

Apply a labelled anti-mouse conjugate at final step (FITC,...).

STORAGE**Pure aliquote**

* at -20°C : until expiry date.

* at + 5°C ± 3°C : 3 weeks.

Diluted aliquote

To be used within 24 hours.

Keep refrigerated (+5°C ± 3°C).

PRESENTATION

1ml vial.

FOR VETERINARY USE ONLY / FOR IN VITRO USE ONLY
SYNBIOTICS EUROPE - 2 rue Alexander Fleming - 69367 LYON Cedex 07 – France
Tel : +33 4.72.76.11.11 - Fax : +33 4.72.76.11.10
www.synbiotics.fr info@synbiotics.fr